

WH-861

One Touch Vortex Shaker

Manual

Madell Technology Corp.

<http://www.madelltech.com>



1、 Technical Data

Rotational speed: 2400r/min

Power: AC110V / 220V


Operation: touching / continuous

Dimension: 130×170x155(mm)

Speed: fixed

Diameter:70mm(Rubber)

2、 Front Panel

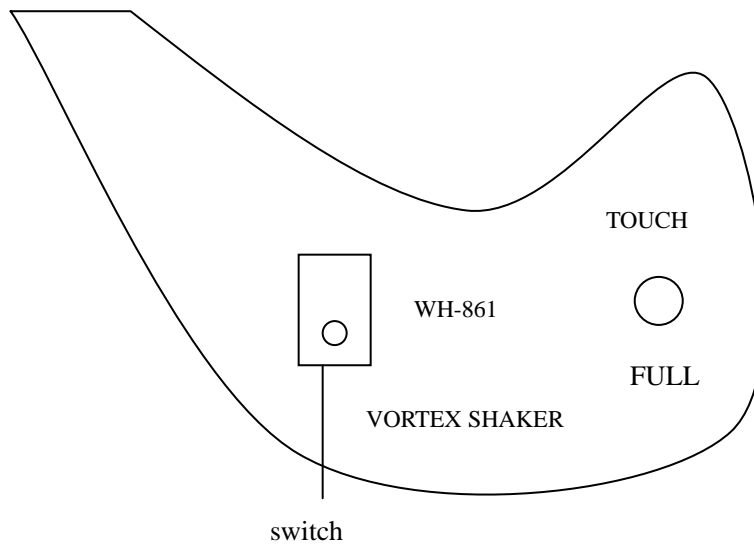

3、 Operations

1. Switch on the power supply.
2. Choose the pattern:
Switch on the "Touch", the mixer is in the touch working condition.
Switch on the "Full", the mixer is in the continuous working condition.
3. Selecting the "Touch" option, the mixer will run when the container is pressed lightly, otherwise it will stop.

Contact:

Madell Technology Corp.
7372 Walnut Ave. Suite V
Buena Park, CA 90620
Toll Free: 1-877-670-9023
<http://www.madelltech.com>