

# ZD-9550 ROCKING SHAKER Manual

Madell Technology Corp.

<http://www.madelltech.com>



# 1. Technical Specifications

Shaking speed: 20-65 rpm

Display: LED

Swinging range:  $\leq 20^\circ$


Platform size: 10.6" × 10.6"

Max load capacity: 10kg

Dimensions: 10.4" × 10.4" × 6"

Supply power: 110V AC, 40W, 50-60HZ

# 2. Front Panel

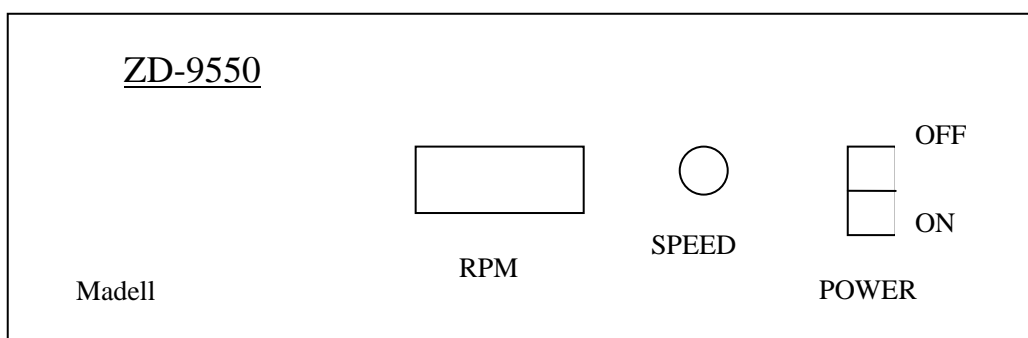

# 3. Operations

1. Connect the power supply; turn on the "Power" switch. The power LED is on to show that the shaker is ready.

2. Adjust the shaking speed: turn the "speed" knob clock-wise makes the shaker faster, counterclock wise makers it slower. The "RPM" display changes accordingly to reflect the new speed.

3. Adjust the swinging angle: this function is located in the middle of the shaker' s back panel. **Before adjusting, turn off the power supply and make sure that the shaker has stopped completely. Otherwise, it may cause serious finger harms.** Loose the roller screw counterclock wise, turn the angle adjustment handle to the proper position, and tighten the screw.

Contact:

Madell Technology Corp.

7372 Walnut Ave. Suite V

Buena Park, CA 90620

USA

Toll Free: 1-877-670-9023

<http://www.madelltech.com>